

ДЕРЖАТЕЛИ 13мм ФИЛЬТРОВ



Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37

Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97

Тверь (4822)63-31-35
Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://millipore.nt-rt.ru> || <mailto:mer@nt-rt.ru>

Swinny Stainless Steel 13 mm Filter Holder


Cat. No. XX3001200

Introduction

The Swinny Stainless Steel 13 mm Filter Holder is a syringe-operated filter holder designed to sterilize or clarify small volumes of liquids by pressure filtration. It can be used to sterilize tissue culture media or clarify HPLC samples, solvents, etc. The filter holder can be autoclaved with the filter in place for clarification/sterilization applications that do not require removing the filter intact for further use.

Filter Loading

1. Unscrew the inlet cover from the outlet base.
2. Remove the O-ring, support screen, and flat gasket.
3. Inspect the gasket, O-ring, and sealing surfaces for damage; imperfections can cause leaks during filtration. Replace damaged parts.
4. Place the flat gasket in the base and place the support screen on top of it.
5. Using smooth-tip forceps, place a 13 mm diameter filter on the support screen, then place the O-ring on top of the filter. If a depth-type prefilter (10 mm diameter) is required to extend the life of the final filter, center the prefilter on top of the final filter before adding the O-ring.
6. Screw the cover down into the base and finger tighten firmly so that the filter is securely sealed.

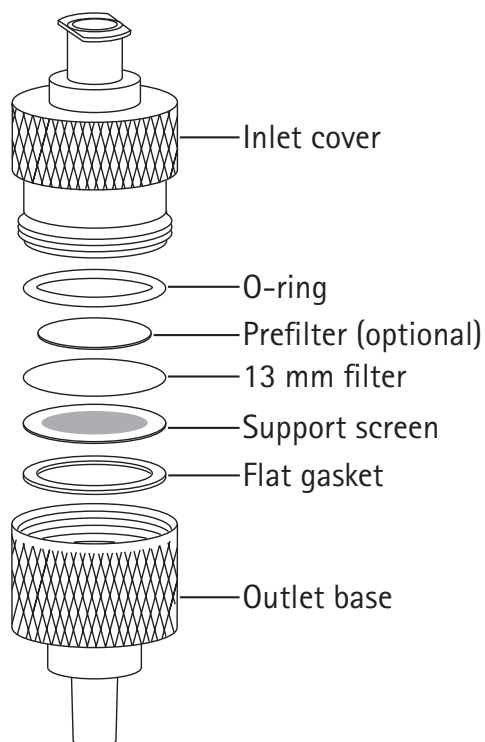

Specifications

Maximum pressure	6.9 bar (100 psi) inlet and differential
Filter diameter	13 mm diameter
Prefilter diameter	10 mm diameter (thick depth prefilter)
Effective filter area	Approximately 0.8 cm ² (0.13 in ²)
Weight	23 g (0.8 oz)
Connections	Female Luer-Lok™ inlet, male Luer-slip outlet
Dimensions	Diameter: 16.0 mm (0.63 in.) Length: 36.6 mm (1.44 in.)
Materials of construction	Inlet cover: 316 stainless steel Outlet base: 316 stainless steel Support screen: 316 stainless steel Flat gasket: polytetrafluoroethylene (PTFE) O-ring: PTFE

Statement Regarding Compliance with the Pressure Equipment Directive, 97/23/EC

EMD Millipore Corporation certifies that this product complies with the European Pressure Equipment Directive, 97/23/EC of 29 May 1997. This product is classified under Article 3 § 3 of the Pressure Equipment Directive. It has been designed and manufactured in accordance with sound engineering practices to ensure safe use. The product is accompanied by user instructions and bears markings to permit identification of EMD Millipore Corporation as the manufacturer or authorized representative of this product within the European Community. In compliance with Article 3 § 3 of the Pressure Equipment Directive, this product does not bear the CE mark.

Ordering Information

This section lists the catalogue numbers for Swinny 13 mm Filter Holder, replacement parts, and accessories.

Description	Catalogue Number
Swinny Stainless Steel 13 mm Filter Holder (includes inlet cover [1], outlet base [1], O-ring [1], support screen [1], flat gasket [1])	XX3001200
Swinny 13 mm Filter Holder Replacement Parts Kit (includes O-ring [4], flat gasket [4], filter holder support screen [4])	XX30012RK
Support screen (1/pk)	XX3001210
Flat gasket, PTFE (10/pk)	XX3001202
Filter forceps, stainless, blunt end (3/pk)	XX6200006P
Plastic syringe, 20 mL capacity (12/pk)	XX1102012
Plastic syringe, 50 mL capacity (5/pk)	XX1105005

Notice

The information in this document is subject to change without notice and should not be construed as a commitment by EMD Millipore Corporation (“Millipore”) or an affiliate. Neither EMD Millipore Corporation nor any of its affiliates assumes responsibility for any errors that may appear in this document.

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37

Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97

Тверь (4822)63-31-35
Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://millipore.nt-rt.ru> || <mailto:mer@nt-rt.ru>