

МИКРОФЛЮИДНАЯ ДРОЖЖЕВАЯ ПЛАТА CELLASIC ONIX Y04D-02



Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37

Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97

Тверь (4822)63-31-35
Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://millipore.nt-rt.ru> || <mailto:mer@nt-rt.ru>

CellASIC® ONIX Y04D-02 Microfluidic Yeast Plate

For research use only. Not for use in diagnostic procedures.

Introduction

The CellASIC® ONIX Y04D-02 Microfluidic Plate is a 4-chamber cell culture plate designed for use with the CellASIC® ONIX2 Microfluidic System and ONIX2 Manifolds for enabling perfusion-based, long-term, live-cell analysis with solution switching. This bio-inspired plate provides a controlled and dynamic microenvironment for cells. The easy-to-use format and superior technology redefine the standard for microfluidics-based experimentation.

Applications

- Time-lapse analysis of yeast cells
- Long-term continuous perfusion experiments
- Solution exchange experiments (induction, inhibition, drug dosing, etc.)
- Comparison up to 4 different cell types or exposure conditions (media components) in parallel
- Cell division tracking (follow mother/daughter cells over generations)
- Temperature and gas atmosphere control (temperature shift, anoxic conditions, etc.)

Plate Description

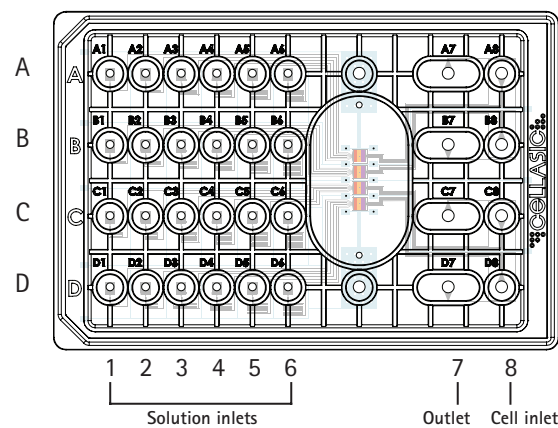


Figure 1. Plate configuration

The Y04D plate has 4 independent units (A–D), each with 6 inlet wells (1–6), a cell inlet (8), and a large outlet well (7). Each row of wells (A–D) addresses the corresponding culture chamber. The plate is shipped prepped with a PBS (phosphate-buffered saline) solution, which can be replaced with a buffer of choice prior to experiment. Each chamber has trap regions of 5.0, 6.0, and 7.0 μm in height to hold cells in a single focal plane during long-term analysis. The plate is for single use only.

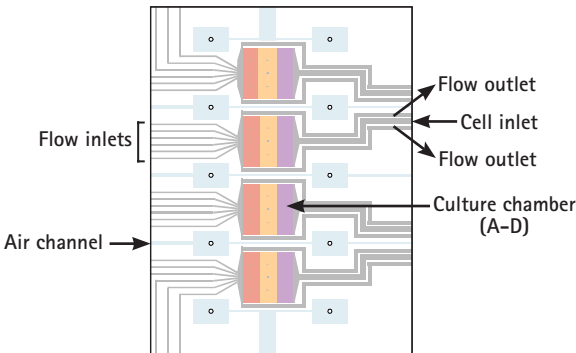


Figure 2. Chamber viewing window

All four culture chambers are located under a single viewing window to minimize travel distance for high-magnification phase objectives.

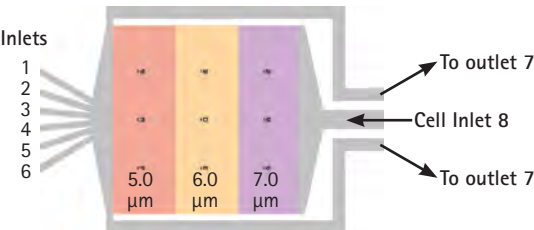


Figure 3. Culture chamber

The culture chamber is 3.0 \times 3.0 mm in area with trap heights of 5.0, 6.0, and 7.0 μm . Nine position markers indicate unit number and relative position. The inlet/outlet functions and minimum/maximum volumes for each culture unit are listed below.

	Function	Minimum Volume (μL)	Maximum Volume (μL)
Inlet 1	Inlet for solution switching	50	350
Inlet 2	Inlet for solution switching	50	350
Inlet 3	Inlet for solution switching	50	350
Inlet 4	Inlet for solution switching	50	350
Inlet 5	Inlet for solution switching	50	350
Inlet 6	Inlet for solution switching	50	350
Inlet 8	Cell inlet for loading cells into culture chamber	50	350
Outlet 7	Accepts flow-through from culture chamber	50	600

Cell Trapping Mechanism

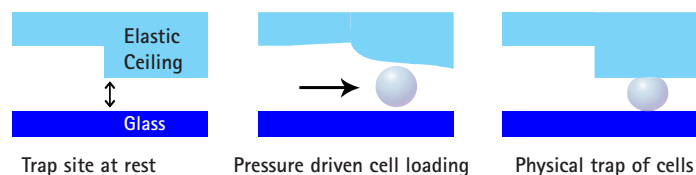


Figure 4. Cell trapping mechanism

The microfabricated chamber gently holds cells against the glass viewing surface to maintain a single focal plane during perfusion analysis experiments. The Y04D plate has trap heights of 5.0, 6.0, and 7.0 μm .

Manifold Description

The CellASIC® ONIX2 heated (CAX2-MXT20) or basic (CAX2-MBC20) manifolds connect the microfluidic plate to the CellASIC® ONIX2 Microfluidic System.

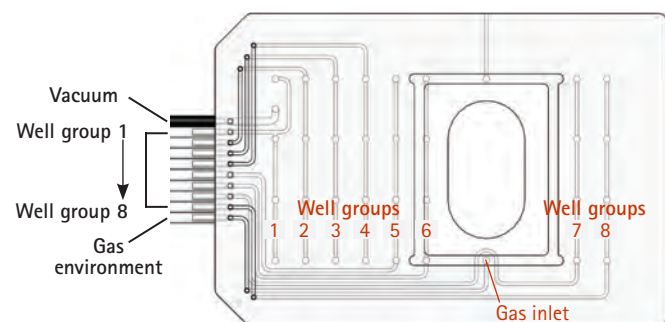


Figure 5. Lines to CellASIC® ONIX2 Microfluidic System

Flow control is achieved using air pressure above the liquid in each well. Multiple wells on a plate are grouped together and addressed by a single pneumatic line via the manifold. Each set of wells is called a "well group". A vacuum line is used to seal the plate to the manifold, and a gas line enables atmospheric control.

Flow Properties

The flow properties of wells 1–6 are shown in Figure 6. The figure shows the flow rate out of the well as a function of pressure. If more than one channel is pressurized, multiply the well flow rate by the number of pressurized channels to derive the overall flow rate.

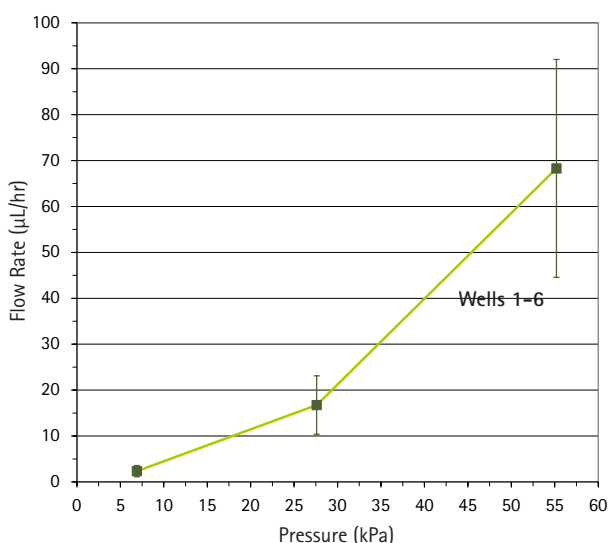


Figure 6. Flow rate for wells 1–6

Plate Storage

Store at room temperature. Do not store in direct sunlight.

Limitations

The plate is incompatible with acetic acid and organic solvents such as acetone, ethanol, and methanol. Plates should be tested for compatibility with other acids or organic solvents prior to use.

Plate Operation

If temperature control is needed, use the CellASIC® ONIX2 Manifold XT (CAX2-MXT20). Refer to the CellASIC® ONIX2 Microfluidic System User Guide for setup instructions.

Plate Preparation

1. Aspirate the PBS solution from wells that will be used for the experiment and add 350 μL of your solution/medium to these wells. Make sure that the unused solution inlet wells are filled with buffer.
2. **Steps 2–4 are optional:** If your experiments require complete removal of PBS, replace the PBS in the solution (1–6) and cell inlet (8) wells with 150 μL of your desired priming solution.
3. Seal the microfluidic plate to the ONIX2 manifold according to the CellASIC® ONIX2 Microfluidic System User Guide.
4. Open the CellASIC® ONIX2 Software, select one of the **New Experiment** options, and find the Y04D plate on the drop down list. On the **Manual Mode** tab (Figure 7), click on the **Run liquid priming sequence** button. Alternatively, on the **Protocol Editor** tab (Figure 8) enter the desired steps and conditions. The recommended pressure and flow time for well groups 1–6 are 34.5 kPa [5 psi] and 5 minutes, respectively. For information on creating a protocol, refer to the CellASIC® ONIX2 Microfluidic System User Guide.

Cell Loading

Pressure-Driven Method Using the CellASIC® ONIX2 Microfluidic System

1. Prepare a yeast/cell suspension of $1.5\text{--}6 \times 10^6$ cells/mL. This concentration may need optimization depending on the yeast strain and desired trapping density.
2. Aspirate solution from cell inlet well 8 and waste outlet well 7.
3. Pipette 50 μL of cell suspension into cell inlet well 8, making sure to cover the hole at the bottom of the well.
4. Seal the microfluidic plate to the ONIX2 manifold according to the CellASIC® ONIX2 Microfluidic System User Guide.
5. Open the CellASIC® ONIX2 Software, select one of the **New Experiment** options, and find the Y04D plate on the drop down list. On the **Manual Mode** tab (Figure 10), click on the **Run cell loading sequence** button. The recommended pressure and flow time for well group 8 are 55.2 kPa [8 psi] and 5 seconds, but you may need to optimize these conditions depending on your cell type/strain and desired trapping density.
6. Assess the loading density on a microscope. If insufficient loading has occurred, repeat the loading protocol.
7. To clear the chamber of untrapped cells, flow one or more inlet well solutions at 34.5 kPa (5 psi) for 5 minutes. On the **Manual Mode** tab, click on the **Run a custom sequence** button or go to **Protocol Editor** to enter the desired parameters. For more information on creating protocols, refer to the CellASIC® ONIX2 Microfluidic System User Guide.
8. Proceed to Cell Culture or Solution Switching sections.

Cell Culture

Cell Culture with CellASIC® ONIX2 Microfluidic System

1. Aspirate solution from wells that will be used for perfusion (wells 1–6). Add 350 μL medium to these wells. Make sure that the unused solution inlet wells are filled with buffer.
2. Seal the microfluidic plate to the ONIX2 manifold according to the CellASIC® ONIX2 Microfluidic System User Guide.

Specifications

Culture Plate Dimensions	
Length x width	127.3 x 85.2 mm (5.0 x 3.4 in.)
Height without lid	14.3 mm (0.6 in.)
Culture Chamber Dimensions	
Length	3.0 mm (0.1 in.)
Width	3.0 mm (0.1 in.)
Trap heights	5.0, 6.0, and 7.0 μm
Glass bottom thickness (#1.5 slide)	170 μm
Plate materials of construction	Polycarbonate, silicone, acrylic, glass

Product Ordering Information

This section lists catalogue numbers for the CellASIC® ONIX products. See Technical Assistance section for contact information.


Description	Catalogue Number	Qty/pk
Microfluidic Plates		
CellASIC® ONIX Plate for Bacteria Cells (4-chamber, trap heights of 0.7, 0.9, 1.1, 1.3, 2.3, and 4.5 µm)	B04A-03-5PK	5
CellASIC® ONIX Gradient Plate for Mammalian Cells (4-chamber)	M04G-02-5PK	5
CellASIC® ONIX Open-top Plate for Mammalian Cells (4-chamber)	M04L-03-5PK	5
CellASIC® ONIX Switching Plate for Mammalian Cells (4-chamber)	M04S-03-5PK	5
CellASIC® ONIX Plate for Haploid Yeast Cells (4-chamber, trap heights of 3.5, 4.0, and 4.5 µm)	Y04C-02-5PK	5
CellASIC® ONIX Plate for Diploid Yeast Cells (4-chamber, trap heights of 5.0, 6.0, and 7.0 µm)	Y04D-02-5PK	5
CellASIC® ONIX Pad Trap Plate (4-chamber, trap height of 4.0 µm)	Y04T-04-5PK	5
CellASIC® ONIX2 Microfluidic System and Manifolds		
CellASIC® ONIX2 Microfluidic System	CAX2-S0000	1
CellASIC® ONIX2 Manifold XT (temperature controlled)	CAX2-MXT20	1
CellASIC® ONIX2 Manifold Basic (no temperature control)	CAX2-MBC20	1
Replacement Parts/Accessories		
CellASIC® ONIX2 Filter Multiconnector (includes filters)	CAX2-AMC00	1
CellASIC® ONIX2 Software USB Drive	CAX2-SSW01	1
CellASIC® ONIX2 Gasket	CAX2-AGK20	1
CellASIC® ONIX2 Self Check Plate	CAX2-ASP20	1
CellASIC® ONIX2 Cleaning Plate	CAX2-ACP20	1
CellASIC® ONIX2 Replacement Filter Pack (9 × 4 mm and 1 × 13 mm Millex® 0.45 µm PTFE filters)	CAX2-AFP00	1
CellASIC® ONIX2 Accessory Fittings (quick-connect gas fitting, 2/pk)	CAX2-ABF00	1
CellASIC® ONIX2 Temperature Calibration Plate	CAX2-ACT20	1

Product Ordering Information, continued

Description	Catalogue Number	Qty/pk
Replacement Parts/Accessories		
CellASIC® ONIX2 Premixed Gas Regulator (for use with 103 L or 112 L gas cylinders with a C10 connection)	CAX2-ABR00	1
CellASIC® ONIX2 Microfluidic Services		
CellASIC® ONIX2 Essential Service Plan	CAX2-ESVC	1
CellASIC® ONIX2 Total Service Plan	CAX2-TSVC	1
CellASIC® ONIX2 Installation	CAX2-INST	1

Notice

The information in this document is subject to change without notice and should not be construed as a commitment by EMD Millipore Corporation ("Millipore") or an affiliate. Neither EMD Millipore Corporation nor any of its affiliates assumes responsibility for any errors that may appear in this document.



Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37

Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97

Тверь (4822)63-31-35
Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://millipore.nt-rt.ru> || <mailto:mer@nt-rt.ru>