

ХРОМАТОГРАФИЧЕСКИЕ КОЛОННЫ ISORAK



Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37

Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97

Тверь (4822)63-31-35
Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://millipore.nt-rt.ru> || <mailto:mer@nt-rt.ru>

IsoPak™ Biochromatography Process Columns

Faster, cleaner, and more reproducible packing and unpacking for process scale chromatography columns

Recognized for its proven chromatographic performance, IsoPak™ columns provide faster, cleaner, and more reproducible packing and unpacking for process scale chromatography columns in purification processes.

The pumped slurry packing and unpacking enables the column to be packed and unpacked without removing the top flow cell. As a result, the operation is contained and avoids exposure of the chromatographic media to the operator and vice versa.

Faster and Reproducible Packing

Columns, even as large as 1400 mm diameter, can typically be packed in 30–60 minutes. The ability to pack a preset bed height at a preset pressure ensures reproducible packing conditions. Given that the slurry pump will automatically stall once the column is fully packed, it cannot over pressurize the bed.

Benefits

- Consistent packing and unpacking provide reproducible results
- Easy to maintain
- Scalable



Scalable

IsoPak™ columns not only provide scalability of column packing, but also through use of our proven flow distribution, scalability of chromatography performance across the range of column diameters. This enables simple and confident scale up of processes from small to large diameter columns. Chromatography performance is equivalent to the best conventional packing methods, enabling processes to be transferred from existing columns with no loss of performance.

Flexible

IsoPak™ columns are easy to pack and are available in a variety of options. These include IsoPak™ valves, antijet, and proven flow cell distribution, providing consistent performance and scalability across a wide range of column diameters.

Availability

IsoPak™ columns are available in a wide range of diameters including 440 mm, 630 mm, 800 mm, 1000 mm, 1200 mm and 1400 mm with column tube in acrylic or stainless steel 316L or 904L. All dimensions are available with variable bed height from 50 to 500 mm or from 500 to 1000 mm.

IsoPak™ Valve

The IsoPak™ three-port valve allows the media slurry to be pumped into the column during packing and the packed bed reslurried in-situ and resuspended media pumped out of the column during unpacking (Figure 1).

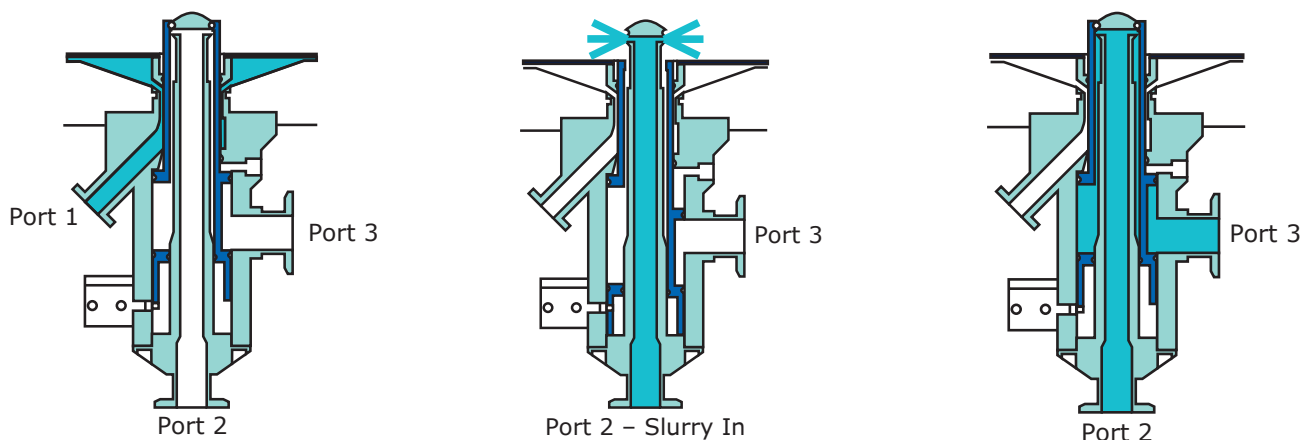


Figure 1: IsoPak™ valve ports operation

Operations Process

In the default or process (run) position, the sleeve is fully extended closing off the slurry ports (Ports 2 and 3) from the column interior.

Packing/Unpacking

When packing, the valve sleeve is retracted, opening Port 2 to the column interior. Media slurry can then be pumped through the open port to pack the column. Port 3 is closed at the slurry transfer skid, preventing slurry exiting the column during packing.

Valve Cleaning

Once packed, the sleeve is moved back to the process position, closing the port. In this position the two slurry ports communicate, enabling the slurry line to be flushed and cleaned in place as required. Once the valve is returned to the process position, the pneumatic lines can be disconnected. Unlike some alternative designs, a key feature of the IsoPak™ valve is that it is not retracted from the bed after packing. This avoids the possibility of voids forming, which may decrease chromatography performance and, in the case of some chromatography media, act as initiation points for bed collapse.

Choose the Best Seal Option for Your Specific Application Needs

Since introducing the Minimum Dead Space (MDS) Inflatable Seal (Air or Liquid Filled), we have demonstrated the efficiency of this sealing design. MDS sealing has higher hygienic performance and its proposed with 2 possible inflating media, compressed air or clean water with 20% ethanol.

Pore Size Bed Support


The bed support, which is available for 10 µm, 20 µm and 30 µm, consists of stainless steel mesh screens that enhance flow distribution efficiency. Unlike plastic support, the screen allows the passage of air without requiring pre-wetting.

Pressure Drop

The graphs below show pressure drop data from IsoPak™ columns fitted with 1 inch IsoPak™ valves. Measurements were made with water at 20 °C. The empty columns were filled with water, tubing was attached to the inlet and outlet process ports, and a flow of water was set up to simulate flow through a

packed bed. The pressure difference between inlet and outlet process ports was recorded. Figure 2 therefore shows the pressure drop due to the column alone. Actual pressure drop during operation will be dependent on the additional pressure drop due to the packed bed.

Pressure drop versus volumetric flowrate

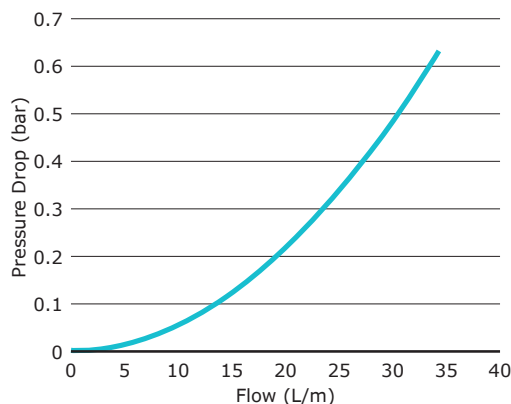

Pressure drop versus linear flowrate

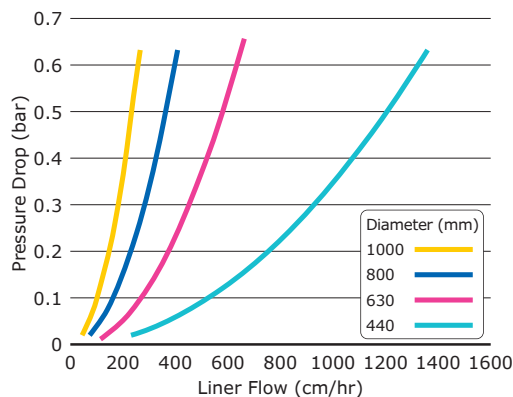


Figure 2: Pressure drop data for IsoPak™ columns fitted with 1 inch IsoPak™ valves

Specifications

IsoPak™ Process Columns

Column Tube Diameter (mm)	440 (441 mm ID), 630 (628 mm ID), 800 (800 mm ID), 1000(990 mm ID), 1200 (1200 mm ID), 1400 (1380 mm ID)
Column Tube Height (mm)	500 or 1000
Maximum Operating Pressure	3.15 bar (45.7 psig)
Operating Temperature (°C)	2 to 30
Material of Construction (Column)	
Column Tube	Acrylic or 316L (1.4404) or 904L (1.4539) stainless steel
Flow Cell	Polypropylene
Bed Support	316L (standard) or 904L stainless steel, 10 µm, 20 µm (standard), 30 µm
Seals	EPDM
Metal Parts	316L or 904L stainless steel (optional); wetted parts are electropolished
Internal Surface Finishing (µm)	
– Column Tube	0.4
– IsoPak™ Valve	0.4
– Tubing	0.4
Regulatory	All wetted polymeric material conforms to FDA 21 CFR 177 and/or have passed USP <88> Biological Reactivity Class VI test and are supplied with relevant material certificates. Pressure Vessel Category: 1
Connection Size	
Process Connection	3/4 in. sanitary clamp for columns diameter ≤1000 mm; 1 in sanitary clamp for column diameter >1000 mm
Slurry Connection	1 in. sanitary clamp for column diameter ≤1000 mm; 1 1/2 in sanitary clamp for column diameter >1000 mm
Dimensions	
Height	For information concerning the dimension and weight of the IsoPak™ column, please consult the arrangement drawings

Antijet column do not have installed Isopak™ valves and are packed manually

Ordering Information

Columns with MDS Air Seal

Product	IsoPak™ Valve	Antijet*	Catalogue No.				Cross Sectional Area (cm²)	Adjustable Height (mm)	Adjustable Capacity (L)	General Arrangement Drawing
			SST 904 L Tube + SST Wetted Parts in 904L	SST 316L Tube + SST Wetted Part in 316L	Acrylic Tube + SST Wetted Parts in 316L	Acrylic Tube + SST Wetted Parts in 904L				
440 x 500 x 450	•		94445321	94443321	94441321	94444321	1520	50–500	8–76	57454WE
		•	94445421	94443421	94441421	94444421				
440 x 1000 x 500	•		94445322	94443322	94441322	94444322	1520	500–1000	76–152	57454WE
		•	94445422	94443422	94441422	94444422				
630 x 500 x 450	•		94635321	94633321	94631321	94634321	3119	50–500	16–156	57459WE
630 x 500 x 450		•	94635421	94633421	94631421	94634421				
630 x 1000 x 500	•		94635322	94633322	94631322	94634322	3119	500–1000	156–312	57459WE
630 x 1000 x 500		•	94635422	94633422	94631422	94634422				
800 x 500 x 450	•		94805321	94803321	94801321	94804321	5027	50–500	25–251	57465WE
800 x 500 x 450		•	94805421	94803421	94801421	94804421				
800 x 1000 x 500	•		94805322	94803322	94801322	94804322	5027	500–1000	251–502	57465WE
800 x 1000 x 500		•	94805422	94803422	94801422	94804422				
1000 x 500 x 450	•		94105321	94103321	94101321	94104321	7697	50–500	39–393	57469WE
1000 x 500 x 450		•	94105421	94103421	94101421	94104421				
1000 x 1000 x 500	•		94105322	94103322	94101322	94104322	7697	500–1000	393–785	57469WE
1000 x 1000 x 500		•	94105422	94103422	94101422	94104422				

Columns with MDS Liquid Seal

Product	IsoPak™ Valve	Antijet	Catalogue No.				Cross Sectional Area (cm²)	Adjustable Height (mm)	Adjustable Capacity (L)	General Arrangement Drawing
			SST 904 L Tube + SST Wetted Parts in 904L	SST 316L Tube + SST Wetted Part in 316L	Acrylic Tube + SST Wetted Parts in 316L	Acrylic Tube + SST Wetted Parts in 904L				
440 x 500 x 450	•		94445331	94443331	94441331	94444331	1520	50–500	8–76	57454WE
440 x 500 x 450		•	94445431	94443431	94441431	94444431				
440 x 1000 x 500	•		94445332	94443332	94441332	94444332	1520	500–1000	76–152	57454WE
440 x 1000 x 500		•	94445432	94443432	94441432	94444432				
630 x 500 x 450	•		94635331	94633331	94631331	94634331	3119	50–500	16–156	57459WE
630 x 500 x 450		•	94635431	94633431	94631431	94634431				
630 x 1000 x 500	•		94635432	94633332	94631332	94634332	3119	500–1000	156–312	57459WE
630 x 1000 x 500		•	94635432	94633432	94631432	94634432				
800 x 500 x 450	•		94805331	94803331	94801331	94804331	5027	50–500	25–251	57465WE
800 x 500 x 450		•	94805431	94803431	94801431	94804431				
800 x 1000 x 500	•		94805332	94803332	94801332	94804332	5027	500–1000	251–502	57465WE
800 x 1000 x 500		•	94805432	94803432	94801432	94804432				
1000 x 500 x 450	•		94105331	94103331	94101331	94104331	7697	50–500	39–393	57469WE
1000 x 500 x 450		•	94105431	94103431	94101431	94104431				
1000 x 1000 x 500	•		94105332	94103332	94101332	94104332	7697	500–1000	393–785	57469WE
1000 x 1000 x 500		•	94105432	94103432	94101432	94104432				
1200 x 500 x 450	•		94125331	94123331	94121331	94124331	11308	50–500	57–565	57475WE
1200 x 500 x 450		•	94125431	94123431	94121431	94124431				
1200 x 1000 x 500	•		94125332	94123332	94121332	94124332	11308	500–1000	565–1130	57475WE
1200 x 1000 x 500		•	94125432	94123432	94121432	94124432				

Columns with MDS Liquid Seal

Product	IsoPak™ Valve	Antijet	Catalogue No.				Cross Sectional Area (cm²)	Adjustable Height (mm)	Adjustable Capacity (L)	General Arrangement Drawing
			SST 904 L Tube + SST Wetted Parts in 904L	SST 316L Tube + SST Wetted Part in 316L	Acrylic Tube + SST Wetted Parts in 316L	Acrylic Tube + SST Wetted Parts in 904L				
440 x 500 x 450	•		94445331	94443331	94441331	94444331	1520	50–500	8–76	57454WE
440 x 500 x 450		•	94445431	94443431	94441431	94444431				
440 x 1000 x 500	•		94445332	94443332	94441332	94444332	1520	500–1000	76–152	57454WE
440 x 1000 x 500		•	94445432	94443432	94441432	94444432				
630 x 500 x 450	•		94635331	94633331	94631331	94634331	3119	50–500	16–156	57459WE
630 x 500 x 450		•	94635431	94633431	94631431	94634431				
630 x 1000 x 500	•		94635432	94633332	94631332	94634332	3119	500–1000	156–312	57459WE
630 x 1000 x 500		•	94635432	94633432	94631432	94634432				
800 x 500 x 450	•		94805331	94803331	94801331	94804331	5027	50–500	25–251	57465WE
800 x 500 x 450		•	94805431	94803431	94801431	94804431				
800 x 1000 x 500	•		94805332	94803332	94801332	94804332	5027	500–1000	251–502	57465WE
800 x 1000 x 500		•	94805432	94803432	94801432	94804432				
1000 x 500 x 450	•		94105331	94103331	94101331	94104331	7697	50–500	39–393	57469WE
1000 x 500 x 450		•	94105431	94103431	94101431	94104431				
1000 x 1000 x 500	•		94105332	94103332	94101332	94104332	7697	500–1000	393–785	57469WE
1000 x 1000 x 500		•	94105432	94103432	94101432	94104432				
1200 x 500 x 450	•		94125331	94123331	94121331	94124331	11308	50–500	57–565	57475WE
1200 x 500 x 450		•	94125431	94123431	94121431	94124431				
1200 x 1000 x 500	•		94125332	94123332	94121332	94124332	11308	500–1000	565–1130	57475WE
1200 x 1000 x 500		•	94125432	94123432	94121432	94124432				
1400 x 500 x 450	•		94145331	94143331	94141331	94144331	14955	50–500	75–748	57476WE
1400 x 500 x 450		•	94145431	94143431	94141431	94144431				
1400 x 1000 x 500	•		94145332	94143332	94141332	94144332	14955	500–1000	748–1495	57476WE
1400 x 1000 x 500		•	94145432	94143432	94141432	94144432				

Ordering Information

IsoPak™ Process Columns

Catalogue No. Structure

94

IsoPak™ Range

Column Diameter

44 = 440 mm
63 = 630 mm
80 = 800 mm
10 = 1000 mm
12 = 1200 mm
14 = 1400 mm

Tube Material

1 = Acrylic tube + SST
wetted parts in 316L
3 = SST 316L Tube + SST
wetted parts in 316L
4 = Acrylic Tube + SST
wetted part in 904L
5 = SST 904L tube + SST
Wetted parts in 904L

Valve or Antijet

3 = IsoPak™ Valve
4 = IsoPak™ Antijet

Adjuster Seal Type

2 = Air filled MDS
3 = Liquid filled MDS

Tube Height

1 = 500 mm
2 = 1000 mm

Spare Parts and Accessories for IsoPak™ Columns

For additional information, please contact your local Technical Service representative.

Slurry Transfer Skids*

Description	Inlet Piping Dimension	Connection	Weight	Dimensions	Catalogue No.
STS-I					
Slurry Transfer Skid 1	OD 1 in. — Dia 25.4 x 1.65	TC 1½ in.	170 kg	L900 x W700 x H1330	MSP88440001
Flexible Hose	OD 1 in. — Dia 25.4 x 1.65	TC 1½ in.			FTP97448*
Pressure gauge + T, STS 1, PRESSURE GAUGE WIKA 233.50/990.22					FTP97450*
Umbilicals UMBILICAL PNEU/ELECT FF+MM 3M LONG					FTP97449*
STS-II					
Slurry Transfer Skid 2	OD 1½ in. — Dia 38.1 x 1.65	TC 1½ in.	190 kg	L900 x W700 x H1330	MSP88440002
Flexible Hose	OD 1½ in. — Dia 38.1 x 1.65	TC 1½ in.			FTP97451*
Pressure gauge + T, STS 2, PRESSURE GAUGE WIKA 233.50/990.22					FTP97452*
Umbilicals UMBILICAL PNEU/ELECT FF+MM 3M LONG					FTP97449*

* To be ordered separately. Required to run the STS



Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37

Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97

Тверь (4822)63-31-35
Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://millipore.nt-rt.ru> || <mailto:mer@nt-rt.ru>