

ВЫСОКОЭФФЕКТИВНЫЕ ФИЛЬТРЫ POLYGARD CN



Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37

Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97

Тверь (4822)63-31-35
Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://millipore.nt-rt.ru> || <mailto:mer@nt-rt.ru>

Polygard® CN Filters

Versatile pleated filters for the clarification and prefiltration of process fluids

Polygard®-CN products are nominally rated filters designed for particle removal applications in liquids and gases. The pleated structure of Polygard®-CN filter media minimizes pressure differentials during the filtration process, and the all-polypropylene construction of these filters conveys low levels of extractables and broad chemical compatibility. Downstream filter performance is enhanced due to efficient particle and microorganism reduction.

Benefits

- Nominal particle retention rating
- Low pressure drop
- Low extractables levels
- Broad range of micron ratings to match an array of particle removal applications
- Designed for rigorous process conditions and broad chemical compatibility
- Ideal for designing scalable solutions from bench top to full-scale manufacturing



Regulatory Compliance

Polygard®-CN filters are designed, developed, and manufactured in accordance with a Quality Management System approved by an accredited registering body to an ISO® 9000 Quality Systems Standard and are shipped with a Certificate of Quality. Each Opticap® XL capsule and cartridge filter is supported by a Validation Guide for compliance with regulatory requirements. For traceability and easy identification, each filter is marked with identifying characteristics.

Multiple Formats Available

Polygard®-CN filters are available in three formats, seven pore sizes, and multiple configurations that vary by filtration area and the type of inlet/outlet connection.

Media Types

Polygard®-CN (nominal)

- 0.3 µm
- 0.6 µm
- 1.2 µm
- 2.5 µm
- 5.0 µm
- 10.0 µm
- 30.0 µm

Filter Formats

- OptiScale® small-scale disposable capsule filters
- Opticap® XL disposable capsule filters
- Cartridge filters

OptiScale® Process Development Screening Tool

OptiScale® disposable capsule filters with Polygard®-CN media provide a convenient small-volume option for process screening and scaling. These "drop in" filters are ideal for evaluating biopharmaceuticals. OptiScale® capsule filters offer speed-to-market strategies for efficiently developing compounds and biotherapeutics.

The OptiScale® disposable capsule is ideally suited for process development and screening. OptiScale® capsules are faster and easier to set up than conventional 47 mm discs.



OptiScale® Filters

Opticap® XL Disposable Capsule Filters

Opticap® XL disposable capsule filters with Polygard®-CN media are available in two filter sizes.

The patented Opticap® XL capsule design allows unparalleled thermal and hydraulic stress resistance in a disposable filter, resulting in reliability, high confidence in the sterility process and improved cleanliness. The unique capsule design with pleated Polygard®-CN media minimizes hold-up volume and reduces production losses.

Convenient and Easy to Use

Opticap® XL capsule filters eliminate the time and the expense associated with assembling, cleaning, and validating stainless steel housings. Adjustable, easy-to-turn, upstream vents and drain valves with o-ring seals and hose barb connections allow for easy process control. Other ease-of-use features include flow directional arrows and ribbed edges for easy gripping even with gloved hands.



Opticap® XL Filters

The Right Size

A range of filter sizes is available to fit all of your application needs, and to allow easy scale-up of your small volume filtration steps to larger, full-scale filtration processes.

The Right Connections

Self-contained and disposable, Opticap® XL capsule filters are supplied with a choice of inlet and outlet connections to optimize your filtration process, including sanitary flanges which enables a higher flow rate, fractional sanitary flanges, and hose barbs.

Cartridge Filters

Polygard®-CN cartridge filters provide high throughput and minimal differential pressure. Cartridges are robust, strong, resilient and are designed to withstand multiple steam-in-place cycles.

A full range of filter sizes is available to suit your application requirements. A variety of connection options are offered for easy adaptation to existing housings.




Cartridge Filters

Specifications


OptiScale®				
Nominal Dimensions				
Inlet/outlet:	Flange/hose barb	Flange/flange	Hose barb/hose barb	
Maximum length:	82 mm (3.24 in.)	74 mm (2.91 in.)	94 mm (3.70 in.)	
Diameter:	69 mm (2.75 in.)	69 mm (2.75 in.)	69 mm (2.75 in.)	
Weight:	2.3 oz (67 g)	2.3 oz (67 g)	2.3 oz (67 g)	
Filtration Area	13.8 cm²			
Materials of Construction				
Filter media:	Polypropylene			
Structural components:	Polycarbonate			
Vent cap:	PVDF			
Internal seal rings:	fluoroelastomers			
Housing Vent	Adjustable vent with male luer and female Luer-Lok™ connections on inlet side of device.			
Maximum Inlet Pressure	5.5 bar (80 psi) at 25 °C			
Oxidizable Substances	Capsules meet the requirements of the USP Oxidizable Substance for Sterile Water for Filtration Test after a water flush of ≤ 100 mL.			
Sterilization by Autoclave	May be autoclaved for 3 cycles of 60 minutes at 126 °C.			
Good Manufacturing Practices	These products are manufactured in a facility which adheres to FDA Good Manufacturing Practices.			
Component Material Toxicity	Component materials were tested and meet the criteria of the USP <88> Reactivity Test for Class VI Plastics. This product is non-toxic per the USP <88> Safety Test.			
Indirect Food Additive	All component materials meet the FDA Indirect Food Additive requirements cited in 21 CFR 177-182.			
	Opticap® XL 5	Opticap® XL 10	Cartridge Filters 4-inch	Per 10-inch
Nominal Dimensions				
Maximum length:	21.6 cm (8.5 in.)	33.5 cm (13.2 in.)	—	—
Diameter:	10.7 cm (4.2 in.)	10.7 cm (4.2 in.)	71 mm (2.75 in.)	70 mm (2.75 in.)
Filtration Area	0.18 m² (1.9 ft²)	0.42 m² (4.5 ft²)	0.18 m² (1.9 ft²)	0.42 m² (4.5 ft²)
Materials of Construction				
Filter media:	Polypropylene		Polypropylene	
Supports:	Polypropylene		Polypropylene	
Structural components:	Polypropylene		Polypropylene	
Vent o-rings:	Silicone		—	
O-rings:	—		Ethylene Propylene (Std)	
Vent/Drain	¼ in. hose barb with double o-ring seal		—	
Maximum Inlet Pressure	5.5 bar (80 psi) at 25 °C 2.8 bar (40 psi) at 60 °C 1.0 bar (15 psi) at 80 °C		— — —	
Maximum Operating Temperature	—		80 °C continuous	
Maximum Differential Pressure				
Forward:	4.8 bar (70 psid) at 20 °C			
NVR Gravimetric Extractables	After autoclaving and a 24 hour soak in ASTM® Type 1 reagent grade water at controlled room temperature:			
	≤25 mg	≤55 mg	—	≤50 mg
Bacterial Endotoxin	Aqueous extraction contains < 0.5 EU/mL as determined by the Limulus Amebocyte Lysate (LAL) Test.			
Oxidizable Substances	Capsules meet the requirements of the USP Oxidizable Substances Test after a water flush of:			
	≤2000 mL	≤5000 mL	—	5000 mL
Sterilization by Autoclave	May be autoclaved for 3 cycles of 30 minutes at 126 °C. (Cannot be steam sterilized in-line)		May be autoclaved for 10 cycles of 30 minutes at 126 °C or steam sterilized for 10 cycles for 30 minutes at 126 °C or hot water sanitized at 80 °C maximum for 30 minutes.	
Component Material Toxicity	Component materials were tested and meet the criteria of the USP <88> Reactivity Test for Class VI Plastics. Polygard®-CN filters meet the requirements of the USP <88> Safety Test utilizing a 0.9% sodium chloride extraction.			
Indirect Food Additive	All component materials meet the FDA Indirect Food Additive requirements cited in 21 CFR 177-182.			
European Pressure Equipment Directive	We certify that this product complies with the European Pressure Equipment Directive, 97/23/EC of 29 May 1997. This product has been classified under Article 3 § 3 of the Pressure Vessel Directive. It has been designed and manufactured in accordance with sound engineering practice to ensure safe use. In compliance with Article 3 § 3 of this Pressure Equipment Directive, this product does not bear the CE mark.		—	—

Typical Clean Water Flow Rates


Opticap® XL 5 Capsule with Polygard®-CN Media –
0.3 µm Nominal (KN03)

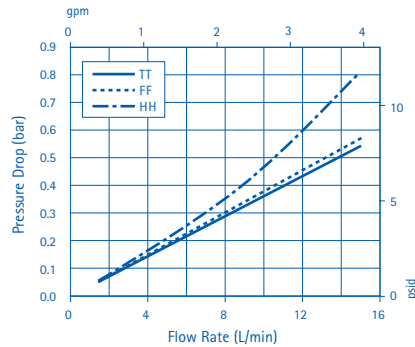


Opticap® XL 5 Capsule with Polygard®-CN Media –
5.0 µm Nominal (KN50)

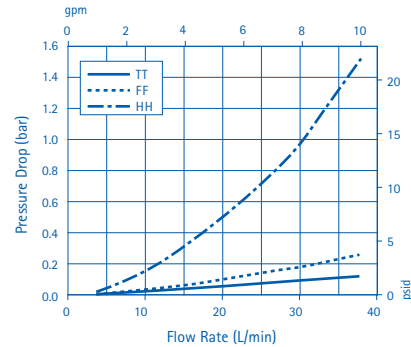

Opticap® XL 10 Capsule with Polygard®-CN Media –
0.3 µm Nominal (KN03)

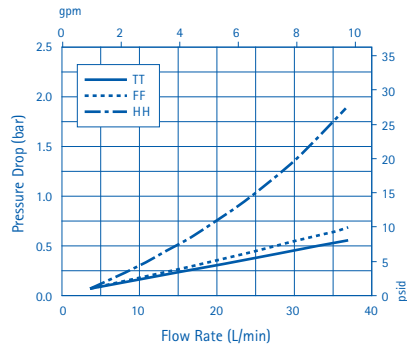

Opticap® XL 10 Capsule with Polygard®-CN Media –
5.0 µm Nominal (KN50)

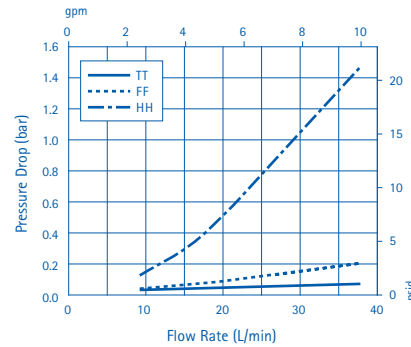

Opticap® XL Legends Refer to Connection Type

TT = 38 mm (1½ in.)
Sanitary Flange Inlet
and Outlet


FF = 19 mm (¾ in.)
Sanitary Flange Inlet
and Outlet

HH = 14 mm (⅝ in.) Hose
Barb Inlet and Outlet

Cartridge Legend Refers to Pore Size


CN03 = 0.3 µm
CN06 = 0.6 µm
CN12 = 1.2 µm
CN25 = 2.5 µm
CN50 = 5.0 µm

Cartridge Filter with Polygard®-CN Media – 10-inch

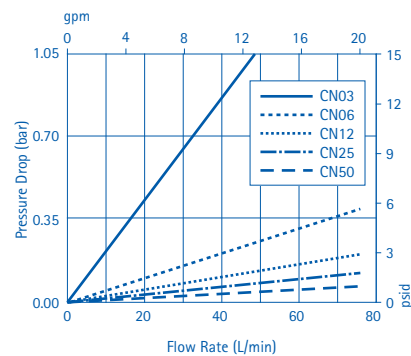


Ordering Information


OptiScale® Capsule Filters

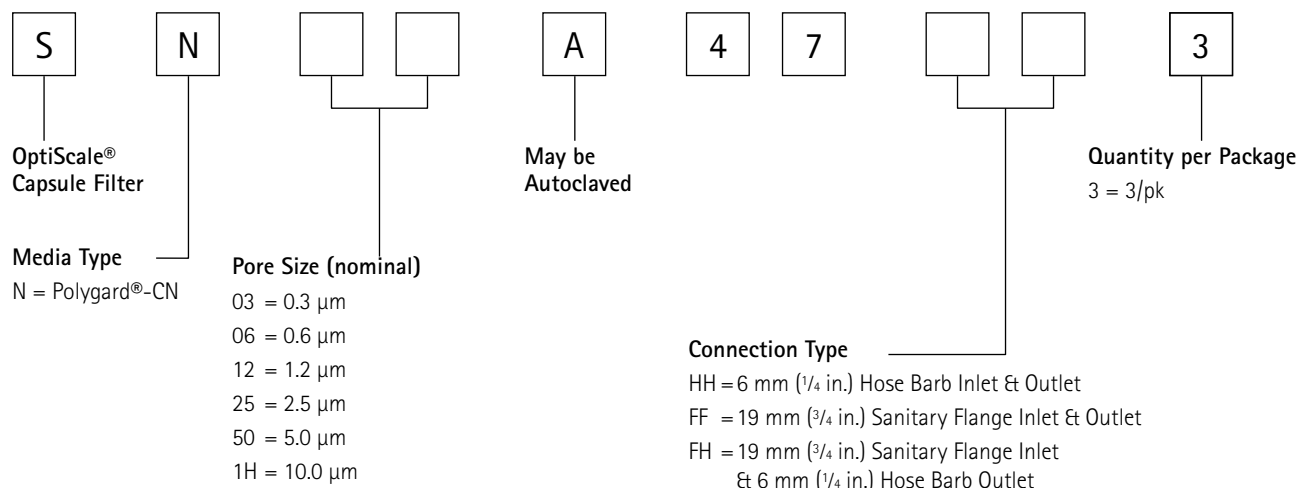

Opticap® XL Capsule Filters

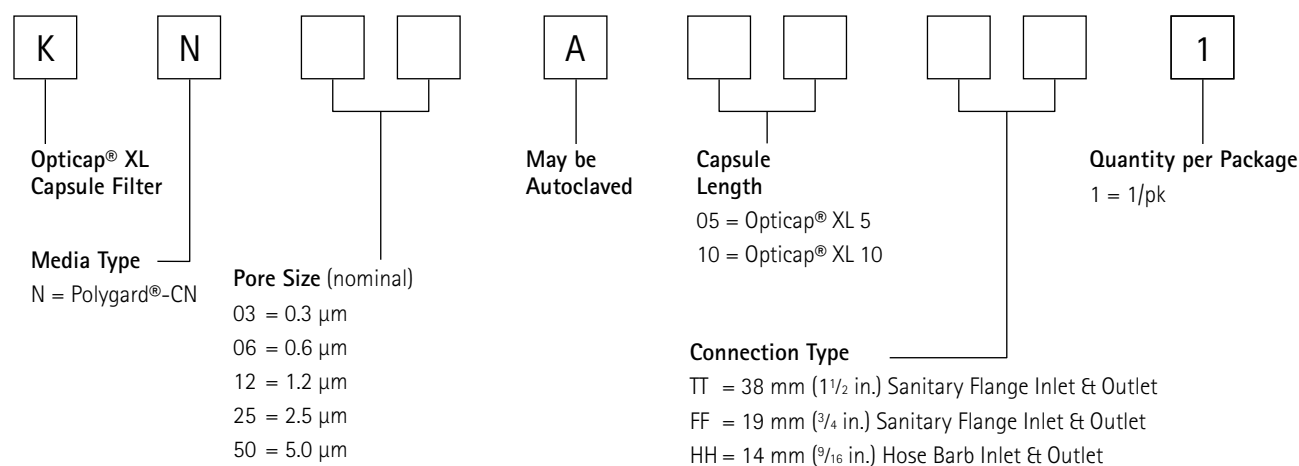

Cartridge Filters

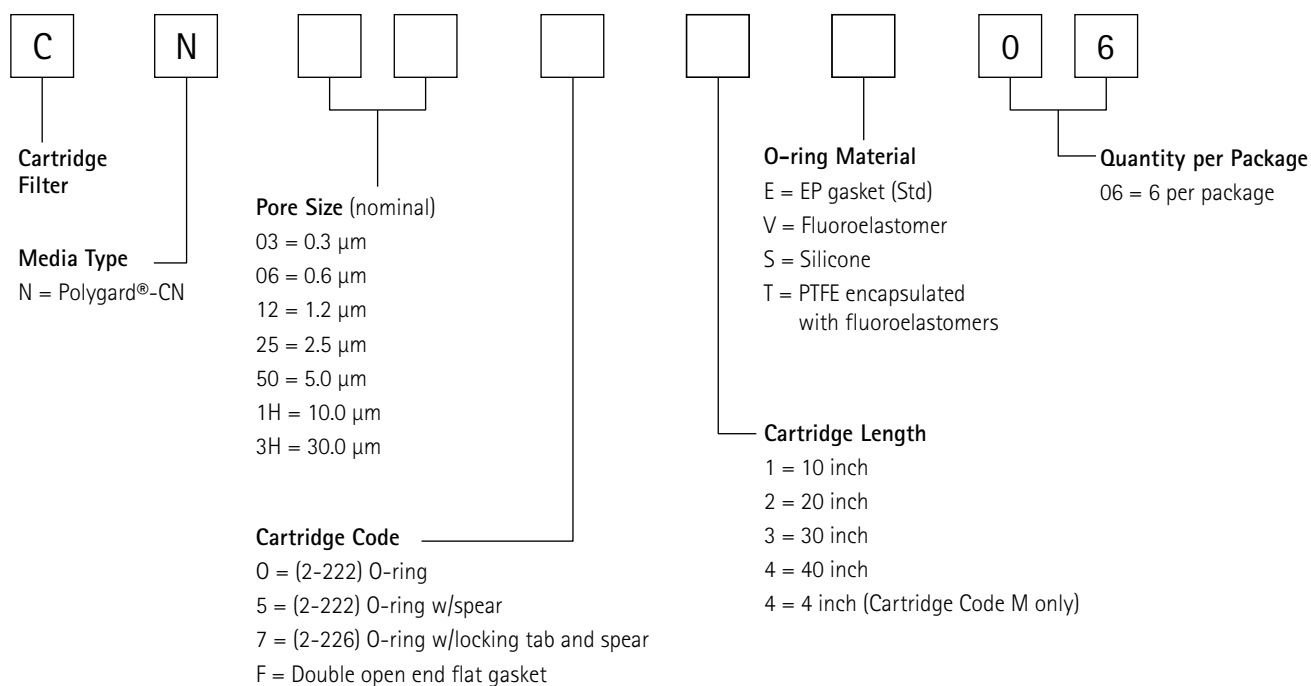



Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37

Пермь (342)205-81-47
Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97

Тверь (4822)63-31-35
Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://millipore.nt-rt.ru> || <mailto:mer@nt-rt.ru>